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removable eye-wash shower with bench support

	cm	Kg	
DE100BR/99	Ø7	2,34	rossa
DE100BV/99	Ø7	2,34	verde
DE100BG/99	Ø7	2,34	gialla

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Sturdy structure obtained from the combination of plastic material joined to an anticorrosive epoxy painted brass body. Soft rubber eyecup with dust cap.

OPERATION: The Goman emergency eye wash showers are made in compliance with the CE, DIN12899, UNI9608 regulations, characterized by a high water flow rate combined with a soft jet that allows immediate washing of the parts of the body accidentally hit by chemicals, acids and corrosive. Indispensable in all high-risk environments and easily identifiable by the red, green or yellow color and the special sign they are equipped with in compliance with the regulations.

INSTALLATION AND OPENING: the installation can be carried out on a bench or on a wall, straight or inclined, single or double, with straight or inclined eyewash. Opening takes place simply by exerting a slight pressure on the operating lever. This lever is equipped with a stop to allow freedom of movement of the hands during emergency operations.

TECHNICAL SPECIFICATIONS CONNECTION: G3/8" to be connected to the water mains. Use only drinking water.

MINIMUM REQUIRED PRESSURE: 2 bar

WATER FLOW: > 6 l/min (as per legislation)

WATER TEMPERATURE: 15÷35°C. Warning: some chemical reactions could be accelerated by an unsuitable temperature. In this case, consult a subject matter expert.

MAINTENANCE: According to the regulations in force in the country where the equipment is used.

REPAIRS: the emergency showers must be repaired only by experienced and qualified personnel. Goman will be relieved of any responsibility for repairs or modifications made by third parties.

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A CONTINUOUS WATER FLOW TO BE OPERATED BY A SINGLE GESTURE ACCORDING TO STANDARD UNI EN 15154-2